1

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The screenshot shows the 'Atelier scientifique' software interface. On the left, the 'Paramétrage' (Parameters) panel is visible, with a 'Temps' (Time) section containing 'Fonction du temps' (Time function) tabs for 'Synchronisation' and 'Mesure'. The 'Mesure' tab is active, showing 'Durée d'acquisition' (Acquisition duration) set to 10 ms, 'Nombre de points' (Number of points) set to 301, and 'Acquisition continue' (Continuous acquisition) unchecked. A green box highlights these settings, with a '2' in a circle. A green dashed arrow points from the 'Mesure' tab to the 'Temps' section, with a '1' in a circle. Below the parameters, there is a 'Sortie analogique S2' section with a green circle '3' around a button. The main window displays a graph with a parabolic curve. A green dashed box highlights the graph area, with a '4' in a circle. A green dashed box highlights the status bar at the bottom, with a '6' in a circle. The status bar shows 'Ordonnées' (Coordinates) and 'Abscisse' (Abscissa). A green dashed box highlights the 'Outils' (Tools) menu, with a '6' in a circle. A green dashed box highlights the 'Fichier' (File) menu, with a '5' in a circle. A green dashed box highlights the 'Graphique' (Graph) menu, with a '5' in a circle. A green dashed box highlights the 'Tableau' (Table) menu, with a '5' in a circle. A green dashed box highlights the 'Compte Rendu' (Report) menu, with a '5' in a circle. A green dashed box highlights the 'Vidéo' (Video) menu, with a '5' in a circle. A green dashed box highlights the 'Afficheur' (Display) menu, with a '5' in a circle. A green dashed box highlights the 'Aide' (Help) menu, with a '5' in a circle.

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